

Synthra MeIplus Research (Catalog No. 003r)

Synthra MeIplus Research is a flexible and completely automated radiosynthesizer for the efficient production of [^{11}C]labeled compounds based on the generation of gas-phase production of [^{11}C]methyl iodide and [^{11}C]methyl triflate. It is specially designed to perform the required multi-step synthesis e. g. for using [^{11}C]propylation. Automating the synthesis is simple, with the easy-to-use configuration software SynthraView. The Synthra MeIplus Research module offers both, fully automatic and manual modes of operation.


Gas Phase Capabilities

- ✓ High specific activities are achieved from in-target produced [^{11}C]CO₂ ranging from 5 Ci/ μmol to 20 Ci/ μmol (Higher specific activities are possible when using methane target).

The [^{11}C]CO₂ produced in target is quantitatively trapped in the stainless steel capillary tubing at -180 °C. Subsequently, the [^{11}C]CO₂ is released into the methane oven where it is converted to [^{11}C]CH₄ by reduction on a Ni-catalyst. The [^{11}C]CH₄ is trapped at -120 °C on Carboxen®. In a successive gas phase reaction the iodination of [^{11}C]CH₄ to [^{11}C]MeI is carried out in a gas phase recirculation system with gaseous I₂ at 730 °C. During circulation [^{11}C]MeI accumulates on a Porapak™ column. Finally, it is released at 200°C and ready for any kind of labeling reaction.

[^{11}C]Labeling Possibilities

- **[^{11}C]Methyl iodide production:** [^{11}C]MeI is ready for release 7 minutes after trapping the [^{11}C]CO₂. The yield for the [^{11}C]methyl iodide formation is under good conditions above 50 % non-decay corrected (ndc).
 - Up to 10 sequential methyl iodide preparations are possible from a single box set-up.
- ✓ **Methyl triflate production:** The [^{11}C]MeI can be converted to [^{11}C]MeOTf by passing through a silver triflate filled column at 180 °C. The conversion yield from methyl iodide is 95 %.
 - Both [^{11}C]MeI and [^{11}C]MeOTf can be used for solid support heterogeneous reactions (e. g. [^{11}C]choline, [^{11}C]methionine) or can be released into the reaction vessel for homogeneous reactions.
- ✓ **Acetate production:** The purified [^{11}C]CO₂ is passed into the reaction vessel for Grignard reactions.

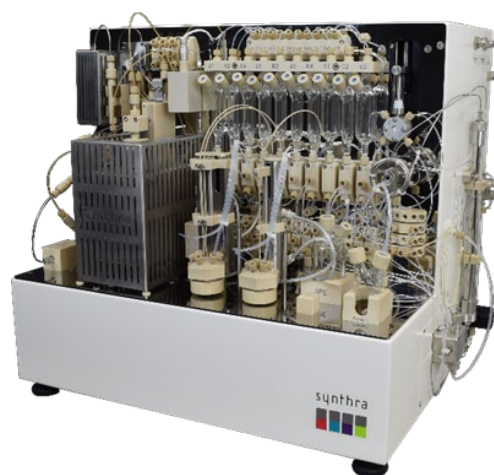

General Features

- ✓ **Heating and cooling capabilities**
 - Nine heating zones
 - Six with cooling capabilities
 - Temperature range: -196 °C – 950 °C
- ✓ **Detectors and controllers**
 - Six shielded radiation detectors
 - Three electronic flow controllers (HCN option: Four flow controllers)
 - Four Pressure sensors
- ✓ **Dispensers and valves**
 - HR-dispenser (up to 50.000 steps, 2.5/5 mL)
 - HPLC pneumatic injection valve (0.5 mL to 2.5 mL sample loop)
 - Five spare valves for customization
 - Chemically inert valves with small dead volume < 35 μL , 5 bar rated
- ✓ **Self-cleaning system option**
- ✓ **Dimensions** (w x d x h): 55 x 50 x 48 cm

Synthra C-11 Family

Product Description and Technical Specifications

synthra



- ✓ **Weight:** approx. 40 kg

Synthesis Features

- ✓ **Two closed reaction vessel** (-196 °C – 200 °C) with integrated cooling to reduce synthesis time
 - 3 mL reaction vessel (minimum volume: 50 µL)
- ✓ **Triflate/column oven** (RT – 200 °C)
- ✓ **Ten reagent vials**
 - Eight small (1 – 3 mL) and two large (10 – 15 mL) volume glass vials for reagents
- ✓ **Three additional cartridge holders**
- ✓ **Built-in preparative radio/UV-HPLC system** for in-process purification and final product collection (max flow: 40 mL/min)
 - Fixed wavelength detector with 255 nm or 280 nm
 - Quaternary gradient
 - One HPLC semi-preparative column
- ✓ **SPE unit** for final product formulation

Additional Synthesis Options

- ➔ **Methane option:** A reduced gas phase suitable for the use of CH₄ target


- ➔ **[¹¹C]CO option** (Catalog No. 003co): After purification, the [¹¹C]CO₂ is released into the column oven for Zn- or Mo-catalyzed reduction to [¹¹C]CO.
- ➔ **[¹¹C]HCN option** (Catalog No. 003hcn): The [¹¹C]CH₄ is released with NH₃ gas into a high temperature area where it undergoes a Pt-catalyzed conversion into [¹¹C]HCN at 950 °C.
- ➔ **Loop option** (Catalog No. 003lo): A heatable and coolable reaction loop is integrated in the synthesis route to reduce synthesis time.
- ➔ **Product solvent evaporator** (Catalog No. 000pse)
- ➔ **Variable wavelength UV detector** (Catalog No. 000vuv)

GMP Features

- ✓ Synthesis files for several [¹¹C]radiotracers available
- ✓ **GMP compliant.** Electronic control and data collection (27/18 channels)
- ✓ **21CFRpart11** & **LIMS** compatible

Terminal Control

- ✓ A laptop (Win 10 Pro) with preinstalled controlling software SynthraView is included
- ✓ Four digital inputs for communication with external devices upon request

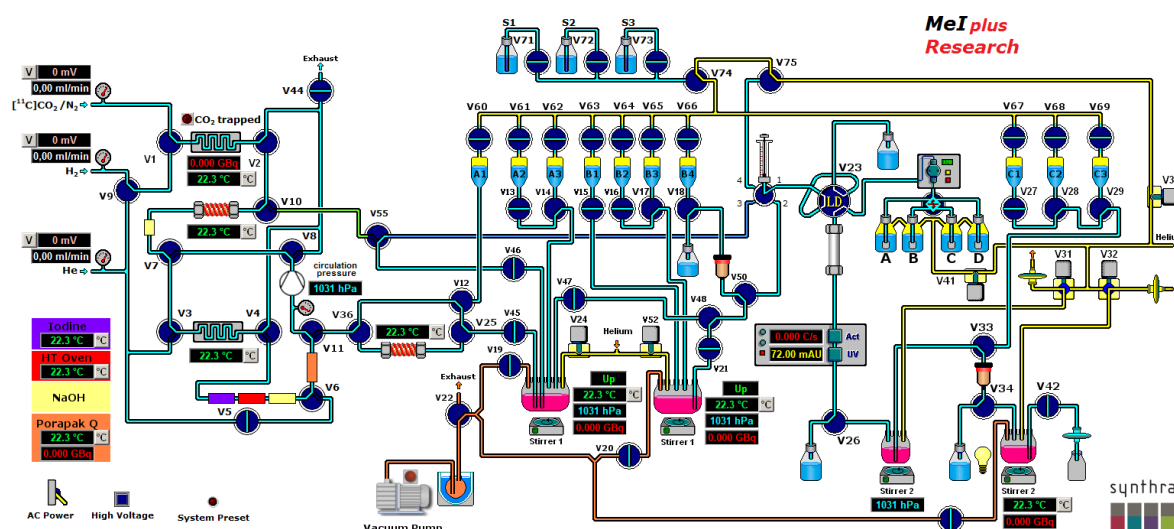

The Graphical User Interface (GUI) of the SynthraView software.

Synthra GmbH reserves the right to modify the information contained herein without prior notice. Copyright © 2022 Printed in the EU